



UNIVERSIDAD
COMPLUTENSE
MADRID

Viral isolation and identification of African swine fever virus in peripheral monocytes

World Organisation for Animal Health (OIE) Reference Laboratory for African swine fever (ASF)

Department of Animal Health, Faculty of Veterinary Medicine, Complutense University of Madrid (UCM).

Veterinary Health Surveillance Center (VISAVET)

Contact:

Prof. José Manuel Sánchez Vizcaíno

jmvizcaino@visavet.ucm.es

Av/ Puerta de Hierro s/n.

28040 Madrid.

Tel: (34) 913944082

Fax: (34) 913943908

Viral isolation and identification of African swine fever virus in peripheral monocytes

1. Virus isolation and identification

1.1 MATERIALS AND REAGENTS

- Donor animal: domestic pig or wild boar (*Sus scrofa*).
- Mechanical defibrinator.
- Adsorbent paper.
- Stopwatch.
- 50 ml tubes.
- CO₂ incubator ($\pm 0.5\%$)/ $37\pm 3^{\circ}\text{C}$.
- Counter chamber [THOMA or NEUBAUER or similar characteristics].
- Eppendorf tubes of 1.5ml.
- Freezer $< -10^{\circ}\text{C}$.
- Freezer $< -70^{\circ}\text{C}$.
- Refrigerator $4\pm 3^{\circ}\text{C}$.
- Glass or plastic pipettes for volumes of 1-25 ml.
- Latex or nitrile gloves.
- Class II laminar flow cabinet.
- 96-well flat-bottom cell culture plate [NUCLON™ Surface, NUNC or similar characteristics].
- Multichannel pipette of 50-300 l.
- Inverted phase contrast cell culture microscope.
- pH meter (0.01 UpH).
- Pipetboy acu or equivalent.
- Incubator with shaker $37\pm 2^{\circ}\text{C}$.
- 250 ml and 500 ml sterile glass bottles.
- Single-channel pipettes 1-10 μl .
- Single-channel pipettes 10-100 μl .
- Single-channel pipettes 10-200 μl .
- Single-channel pipettes 200-1000 μl .

- Tabletop centrifuge with 4°C cooling, Megafuge 1.0R [Heraeus rotor #7570 or similar].
- Vortex.
- Ammonium chloride NH₄Cl [Ref.: 1.01145.1000 (Merck) or similar characteristics]. Store at 4°C.
- Phosphate buffered saline solution (PBS 1x) in tablets [Ref.: 524650-1 (CALBIOCHEM) or similar characteristics]. Store at room temperature.
- Türk's dye. [Ref.: 1.09277.0500 (Merck) or similar characteristics]. Store at room temperature.
- Positive control: ASF virus positive ADH isolate. Store at ≤ -70°C.

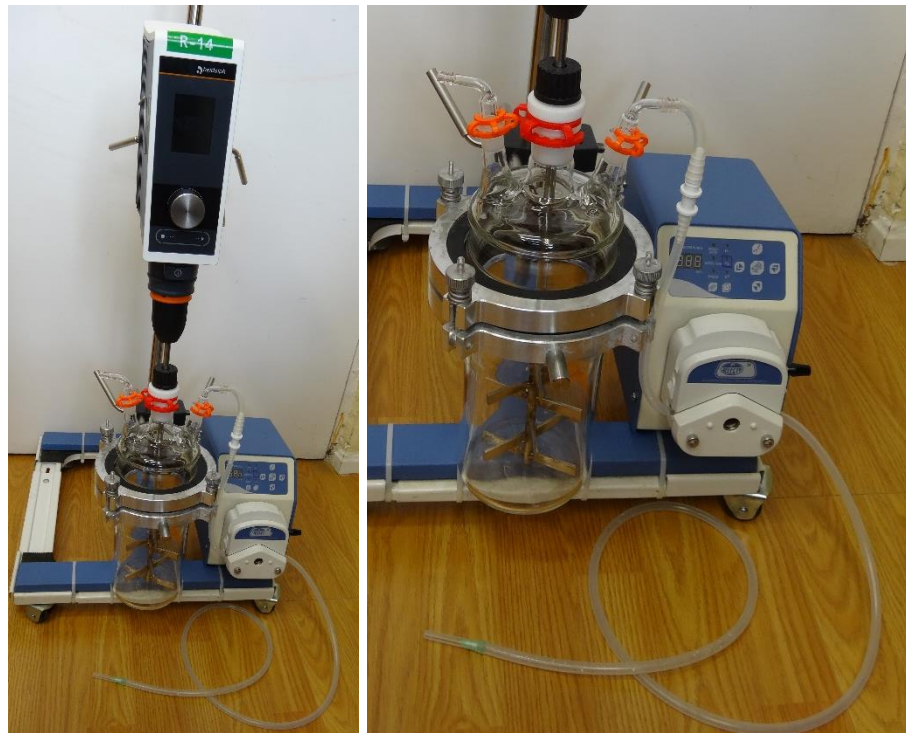


Figure 1. Mechanical defibrinator.

1.2 METHODOLOGY:

1.2.1 EXTRACTION AND CULTURE OF LEUKOCYTES:

- Obtaining defibrinated blood from a donor pig or wild boar:
 - Sedate the animal.
 - Before the blood extraction, introduce the needle with which the extraction will be performed into the tube that connects to the defibrinator, previously started (Figure 1).
 - Perform the blood extraction through the ophthalmic venous sinus.
 - Once the blood collection has been completed, leave the blood inside the defibrillator for 30 minutes with the blades rotating.

- Plate preparation:
 - **Reagents:**
 - Erythrocyte lysis solution: Ammonium chloride NH₄Cl, sterile at 0.83% (8.3gr of NH₄Cl in 1 liter of distilled water).
 - Phosphate buffered saline solution: (PBS 1X, pH ±2): Tablets can be used directly (1 tablet in 1 liter of distilled water) or manufactured as follows:

ClNa- -----	8gr
ClK -----	0.2gr
PO ₄ H ₂ K -----	0.2gr
PO ₄ HNa ₂ -----	1.15gr
Distilled water -----	1 litro
 - Discard the fibrin clot generated by the defibrinator.
 - Aliquot the blood in the defibrinator in 50ml sterile tubes in a laminar flow cabinet.
 - Centrifuge for 30 minutes at 2,500rpm, without brake.
 - We will obtain three phases: serum (culture medium for leukocytes), a thin white layer of leukocytes and a third fraction of erythrocytes.
 - Transfer the serum to a sterile 250 ml jar and use it in the last step as culture medium.
 - Transfer the erythrocytes to a 50 ml tube and store at 4°C.
 - Transfer the leukocytes to a 50 ml tube and add 3 volumes of erythrocyte lysis solution.
 - Incubate on ice for 15 minutes.
 - Centrifuge at 2,000 rpm for 15 minutes, with brake.
 - Discard the supernatant, carefully and add, again, 3 volumes of erythrocyte lysis solution.

- Incubate on ice for 15 minutes.
- Centrifuge at 2,000 rpm for 15 minutes, with brake.
- Discard the supernatant. If the pellet still contains erythrocytes, perform one or two washes with 1X PBS (centrifugations will be at 2,000rpm for 10 minutes).
- Resuspend the leukocyte pellet with 10ml of the previously collected serum.
- Perform the leukocyte count and adjust the leukocyte suspension to a final concentration 10⁷ leukocytes/ml, for this purpose use Turk's solution (5 l leucocyte suspension + 495 µl of Turk's solution).
- Distribute the leukocytes in 96-well plates, 200 µl per well (300,000 cells/well).
- Incubate at 37°C with 5% CO₂ for 3-4 days.

1.2.2 INOCULATION OF SAMPLES AND CONTROLS:

- After 3-4 days, dilute the erythrocytes 1/100 in 1X PBS and inoculate 20 µl of the dilution into each well.
- Perform a 1/10 dilution of each sample in the well to be tested (20 µl/well).
- Inoculate 8 wells per sample, leaving 8 free wells between samples to avoid contamination.
- Inoculate the positive control and negative control.
- Seal the plates with Parafilm.
- Incubate the 96-well plate at 37°C with 5% CO₂ for 5-7 days to check for hemadsorption (HAD) or cytopathic effect (CPE).

1.3 ANALYSIS AND INTERPRETATION OF RESULTS:

- Inoculated wells should be observed daily under the microscope to check for the presence or absence of hemadsorptions (HAD) and/or cytopathic effect.
- The first observation can be made 14-16 hours after inoculation.
- To check for the presence of ADH, plates should be gently shaken before microscopic observation.
- The observation period should be extended to check for the presence of ADH or cytopathic effect up to 7 days after sample incubation.
- After 7 days, viral genome extraction and real-time PCR of the leukocyte culture supernatant from the plate is performed.
- Collection of 10 µl of the sample or biological matrix and a new pass is performed on a new plate with prepared leukocyte culture.

- This procedure is carried out a total of three times, thus obtaining three successive passes of the environmental sample or biological matrix on a leukocyte culture.

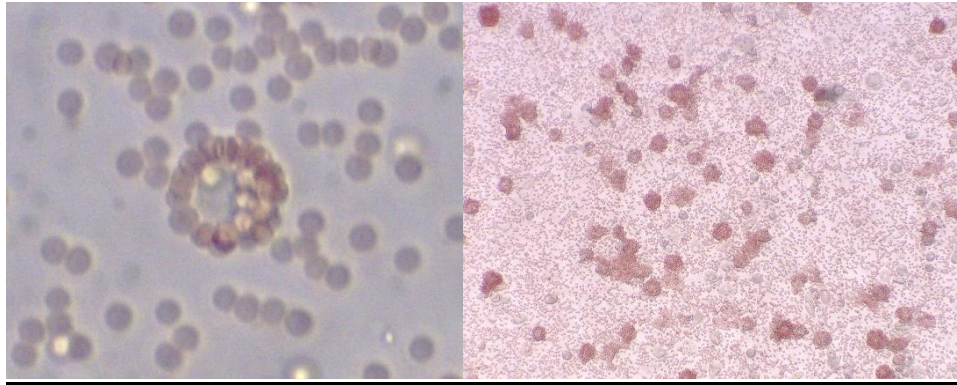


Figure 2. Hemadsorption.

Bibliographic References:

1. Sánchez-Vizcaíno JM, Laddomada A, Arias ML. "African Swine Fever Virus," in *Diseases of Swine* (Hoboken, NJ, USA: John Wiley & Sons, Inc.), 443–452. doi:10.1002/9781119350927.ch25
2. Manual de las pruebas de diagnóstico y de las vacunas para los animales terrestres 2019 de la OIE. Capítulo 2.8.1 OIE, 2019. https://www.oie.int/fileadmin/Home/eng/Health_standards/tahm/3.08.01_ASF.pdf